# SD Series Bench Scales

## User Guide

EN

PT

SV

DA

PL

NL

ES

FR

DE

IT

CZ

HU

# 1. SAFETY INFORMATION

## Definition of Signal Warnings and Symbols

<b>WARNING</b>	For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
<b>CAUTION</b>	For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
<b>Attention</b>	For important information about the product.
<b>Note</b>	For useful information about the product.

## Warning Symbols

General hazard

Electrical shock

## Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

## Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

# 2. INSTALLATION

## 2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.


## 2.2 Connecting Power

### AC Adapter Installation

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

**Attention:** Only use an AC adapter specified by OHAUS.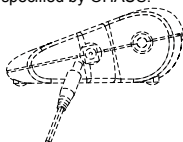


**Battery Installation**

Install the three “C” size (LR14) batteries (not included) with polarity as shown in the battery compartment.

**2.3 Levelling the Instrument**

To level the instrument, adjust the feet/leveling wheel so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.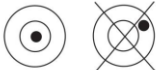

**2.4 Initial Calibration**

For best results, the instrument’s calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

**3. OPERATION**

**3.1 Overview of Display, Home Screen**

**CONTROLS**


**Two Button Keypad**


- **ON/ZERO Off** : A short press turns on the scale when it is off. When the scale is on, a short press zeroes the scale, and a long press turns the scale off.
- **UNITS Cal** : A short press changes the weighing unit and weighing mode.  
A short press initiates countdown during Dynamic Weighing Mode.  
A long press initiates the calibration process.

**Dynamic Weighing**

1. With platform empty, press the **UNITS Cal** button until the indicator • flashes next to the desired measuring unit.
2. Place a mass on the platform. The scale will show “-A-”.  
NOTE: If using a container, press **ON/ZERO Off** to tare the weight and return to zero.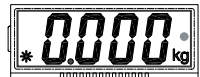

3. With a mass on the platform or container, a short press of UNITS *Cal* will start the averaging countdown from 5 seconds.
4. The averaged weight is then displayed. The indicator • stops flashing when the process is complete.
5. The display will hold until a button is pressed.
6. To repeat the process, press UNITS *Cal* with a mass on the platform or container.
7. To Exit, remove the mass from the platform or container and press UNITS *Cal*.

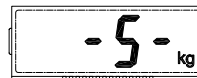## 4. MAINTENANCE

### 4.1 Cleaning

**WARNING:** Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

**Attention:** Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

### 4.2 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website [www.ohaus.com](http://www.ohaus.com) to locate the Ohaus office nearest you.

## 5. TECHNICAL DATA

### Specifications:

Model	SD35	SD75	SD200
Capacity x Readability	35 kg x 0.02 77 lb x 0.05	75 kg x 0.05 165 lb x 0.1	200 kg x 0.1 440 lb x 0.2
Battery power requirements	3 "C" Alkaline batteries (LR14)		
AC Adapter	9-12 Volt DC, 100mA		

Model	SD75L	SD200L
Capacity x Readability	75 kg x 0.05 165 lb x 0.1	200 kg x 0.1 440 lb x 0.2
Battery power requirements	3 "C" Alkaline batteries (LR14)	
AC Adapter	9-12 Volt DC, 100mA	

## 6. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> .

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Ohaus Corporation